


Configuration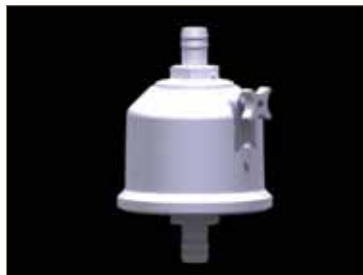

1/2" Hosebarb (Code H)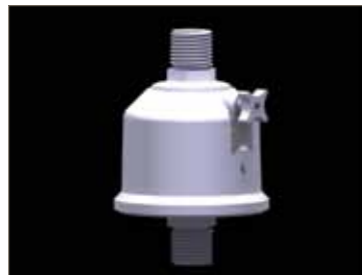

1/4" NPT (Code M)

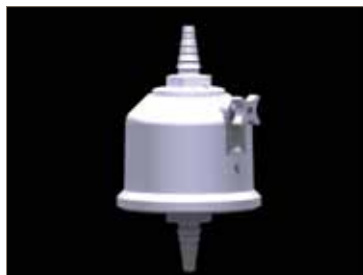
Stepped Hosebarb (Code G)

1" Tri-clamp (Code T)


Gromelle (Code R)

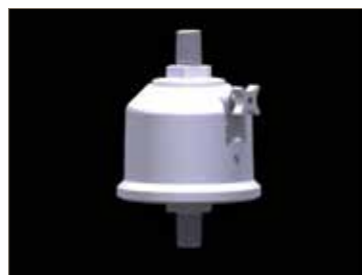
1/2" NPT (Code N)


Walther Male (Code Q)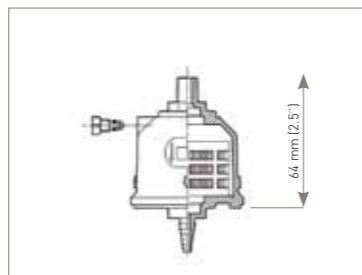

E' Size

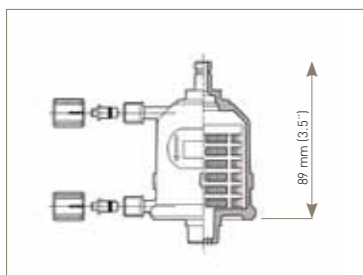B' Size

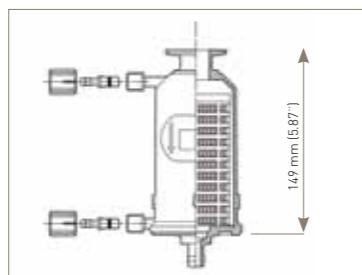A' Size

Sales Offices Worldwide

AE – United Arab Emirates, Dubai
Tel: +971 4 8127100
parker.me@parker.com

AR – Argentina, Buenos Aires
Tel: +54 3327 44 4129

AT – Austria, Wiener Neustadt
Tel: +43 (0)2622 23501-0
parker.austria@parker.com

AT – Austria, Wiener Neustadt
Tel: +43 (0)2622 23501 970
parker.easteurope@parker.com

AU – Australia, Castle Hill
Tel: +61 (0)2-9634 7777

AZ – Azerbaijan, Baku
Tel: +994 50 2233 458
parker.azerbaijan@parker.com

BE/LU – Belgique, Nivelles
Tel: +32 (0)67 280 900
parker.belgium@parker.com

BR – Brazil, Cachoeirinha RS
Tel: +55 51 3470 9144

BY – Belarus, Minsk
Tel: +375 17 209 9399
parker.belarus@parker.com

CA – Canada, Milton, Ontario
Tel: +1 905 693 3000

Tel: +41 (0) 21 821 02 30
parker.switzerland@parker.com

CN – China, Shanghai
Tel: +86 21 5031 2525

CZ – República Checa, Kletcany
Tel: +420 284 083 111
parker.czechrepublic@parker.com

DE – Alemania, Kaarst
Tel: +49 (0)2131 4016 0
parker.germany@parker.com

DK – Denmark, Ballerup
Tel: +45 43 56 04 00
parker.denmark@parker.com

ES – Spain, Madrid
Tel: +34 902 33 00 01
parker.spain@parker.com

FI – Finland, Vantaa
Tel: +358 (0)20 753 2500
parker.finland@parker.com

FR – France, Contamine-sur-Arve
Tel: +33 (0)4 50 25 80 25
parker.france@parker.com

GR – Greece, Athens
Tel: +30 210 933 4450
parker.greece@parker.com

HK – Hong Kong
Tel: +852 2428 8008

HU – Hungary, Budapest
Tel: +36 1 220 4155
parker.hungary@parker.com

IE – Ireland, Dublin
Tel: +353 (0)1 464 6370
parker.ireland@parker.com

IN – India, Mumbai
Tel: +91 22 6513 7081-85

IT – Italy, Corsico (MI)
Tel: +39 02 45 19 21
parker.italy@parker.com

JP – Japan, Tokyo
Tel: +81 3 6408 3901
Tel: +82 2 559 0400

KZ – Kazakhstan, Almaty
Tel: +7 7272 505 800
parker.easteurope@parker.com

LV – Latvia, Riga
Tel: +371 4 745 2601
parker.latvia@parker.com

MX – Mexico, Apodaca
Tel: +52 81 8156 6000

MY – Malaysia, Subang Jaya
Tel: +60 3 5638 1476

NL – The Netherlands, Oldenzaal
Tel: +31 (0)541 585 000
parker.nl@parker.com

NO – Norway, Ski
Tel: +47 64 91 10 00
parker.norway@parker.com

NZ – New Zealand, Mt Wellington
Tel: +64 9 574 1744

PL – Poland, Varsovia
Tel: +48 (0)22 573 24 00
parker.poland@parker.com

PT – Portugal, Leca da Palmeira
Tel: +351 22 999 7360
parker.portugal@parker.com

RO – Romania, Bucharest
Tel: +40 21 252 1382
parker.romania@parker.com

RU – Russia, Moscow
Tel: +7 495 645-2156
parker.russia@parker.com

SA – Republic of South Africa, Kempton Park
Tel: +27 (0)11 961 0700
parker.southafrica@parker.com

SE – Sweden, Spånga
Tel: +46 (0)8 59 79 50 00
parker.sweden@parker.com

SG – Singapore
Tel: +65 6887 6300

SK – Slovakia, Banská Bystrica
Tel: +421 484 162 252
parker.slovakia@parker.com

SL – Slovenia, Novo Mesto
Tel: +386 7 337 6650
parker.slovenia@parker.com

TH – Thailand, Bangkok
Tel: +662 717 8140

TR – Turkey, Istanbul
Tel: +90 216 4997081
parker.turkey@parker.com

TW – Taiwan, Taipei
Tel: +886 2 2298 8987

UA – Ukraine, Kiev
Tel: +380 44 494 2731
parker.ukraine@parker.com

UK – England, Warwick
Tel: +44 (0)1926 317 878
parker.uk@parker.com

US – USA, Cleveland
Tel: +1 216 896 3000

VE – Venezuela, Caracas
Tel: +58 212 238 5422

© 2011 Parker Hannifin Corporation
Stock N°: 17 950 1234
11/11 Rev. 2C

Parker Hannifin Manufacturing Ltd
domnick hunter
Process Filtration - Europe
Durham Road
Birtley, Co. Durham
DH3 2SF, England
phone +44 (0)191 4105121
fax +44 (0)191 4105312
email: dhprocess@parker.com
www.parker.com/processfiltration

Parker Hannifin Corporation
domnick hunter
Process Filtration - North America
2340 Eastman Avenue
Oxnard, California, USA 93030
toll free: +1 877 784 2234
phone: +1 805 604 3400
fax: +1 805 604 3401
email: dhpsales.na@parker.com
www.parker.com/processfiltration

Installation and Operating Instructions for DEMICAP Capsules

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English

Introduction

These guidelines give the correct methods for installing and operating Parker domnick hunter DEMICAP Capsules. If you have any queries please contact our process filtration specialists who will be pleased to discuss your particular filtration requirements or answer any questions you may have. We may also be contacted at any of the addresses given on the reverse of this document or through our worldwide network of distributors.

1. Storage

Store Capsules in a clean and dry environment and avoid placing heavy objects on top of the Capsule box. Capsules should not be stored at temperatures below 5 °C (41 °F) or above 30 °C (86 °F) or in direct sunlight.

1.1 The shelf-life for DEMICAP Capsules is as follows:

Non-gamma irradiated	3 years
Gamma irradiated	2 years
PROPOR PES DEMICAP Capsules	Consult Certificate of Conformance

2. Operating Conditions

Liquid:	Maximum 5 barg (72 psig) @ 40°C
Gas:	Maximum 4 barg (58 psig) @ 40°C

If DEMICAP Capsules are used outside these operating conditions or with fluids that are not compatible with the materials of construction then the product body may fail and result in injury to the operator or damage to equipment.

3. Installation and Operation

Please ensure that the Capsule you have chosen is compatible with the fluid and system before installation and check that the product and part number etched on the base of the Capsule are as required.

NB: If the capsule is to be used in a sterile environment please ensure that the capsule is either sterile packed by Parker domnick hunter or autoclaved prior to use. DEMICAP Capsules can not be steam sterilised in-situ.

A) Liquid Systems

- Follow the arrows on the capsule label and ensure that the Capsule is installed in the correct direction and using the appropriate connections.
- Before passing liquid through the capsule ensure that the vent valves are closed and the Capsule is mounted in a secure position. It is important that the Capsule is full of liquid during operation, so ensure that it is mounted so that the air can be vented in a safe manner.
- The Capsule can be vented by turning the vent nut in an anti-clockwise direction. The valve is retained by the nut so do not undo the nut by more than 1 or 2 turns as this may result in wastage of the liquid within the Capsule or injury to the operator. After long term use of the Capsule, build up of residue on the seats of the vents valves may cause them to stick. Pulling gently on the hosebarbs after undoing the nut by a quarter turn will release them.
NB: If hazardous fluids are to be passed through the Capsule, the vent valves should be connected to an appropriate drain line.
- If the Capsule is connected into a pressurised system make sure that transient pressures do not exceed the maximum operating conditions.

B) Gas Systems

- Follow the arrows on the capsule label and ensure that the capsule is installed in the correct direction and using the appropriate connections. For vent applications flow can be in either direction.
- Gas filters are hydrophobic and may not allow the passage of water. It is also possible that condensation can collect in the Capsule and prevent the passage of gas. The Capsule should be mounted so that any water or condensate can be drained from the vent valves. In some cases it may be necessary to remove the Capsule and dry the membrane by passing dry air or nitrogen through it.

4. Integrity Testing

Some DEMICAP Capsules can be integrity tested either off-line or in-situ. Please contact Parker domnick hunter or its representative for further information on the most appropriate method and equipment. Integrity test parameters are available in the product datasheets.

5. Sterilisation

DEMICAP Capsules can be autoclaved up to the maximum temperature marked on the product. Please contact Parker domnick hunter with the details of your sterilisation regime so we can advise the maximum number of cycles.

- NB:
- The vent / drain valves should be left open during the autoclave cycle
 - DO NOT STEAM STERILISE DEMICAP CAPSULES IN-LINE !

6. Disposal

All Capsules should be disposed of in a safe manner and in line with Health and Safety Guidelines.

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